

Using the autoclave (NL/ENG)

DE AUTOCLAAF GEBRUIKEN

- Haal voor gebruik de binnenpan uit de autoclaaf en beoordeel de hoeveelheid water, dat moet ong. 3,5 liter zijn, of tot de onderkant van de pootjes waarop de binnenpan rust.
- Zet de te steriliseren objecten in de binnenpan. Draai de doppen van flessen losjes op de hals en max 75% vol (om lucht te kunnen laten ontsnappen). Haal evt. parafilm tape eraf voordat je steriliseert.
- Zet de binnenpan terug en sluit het deksel gelijkmatig met de vleugelmoeren.
- Stel de gewenste sterilisatietijd in, wacht 5 seconden en schakel de autoclaaf uit en weer in om de nieuwe tijd te bevestigen.
- De cyclus loopt nu. Als de watertemperatuur 100 graden is klinkt een serie piepjes. Het sterilisatieproces start, de tijd loopt af.
- Nadat de tijd om is klinkt een alarm. Schakel de pan uit. Als de druk gedaald is, open je de pan. **PAS OP, ALLES IS HEET!**
- Na gebruik is het belangrijk om de binnenpan en de inleg helemaal droog te maken, anders gaan deze onderdelen roesten.

USING THE AUTOCLAVE

- Before use, remove the inner pan from the autoclave and assess the amount of water, ideally about 3.5 litres, or to the bottom of the legs on which the inner pan rests.
- Place the objects to be sterilized in the inner pan. Tighten the caps of bottles loosely on the neck and max 75% full (to allow air to escape). Remove the paraffin film before placing Petri dishes in the autoclave.
- Replace the inner pot and close the lid evenly with the wing nuts.
- Set the desired sterilization time, wait 5 seconds and switch the autoclave off and on again to confirm the new time.
- The cycle is now running. When the water temperature reaches 100 degrees, a series of beeps will sound. The sterilization process starts, the time counts down.
- After the time is up, an alarm will sound. Turn off the pan. When the pressure has dropped, open the pan. **BEWARE, EVERYTHING IS HOT!**
- After use, it is important to dry the inner pan and the insert completely, otherwise these parts will rust.

Revision #3

Created 2024-07-15 16:48:32 UTC by Shirley

Updated 2024-07-15 16:57:36 UTC by Shirley